



memmert
Experts in Thermostatics

wiring diagram




compressor-cooled incubator

designation: ICP eco


size: 260

nominal voltage: 230V


drawing number: D39571

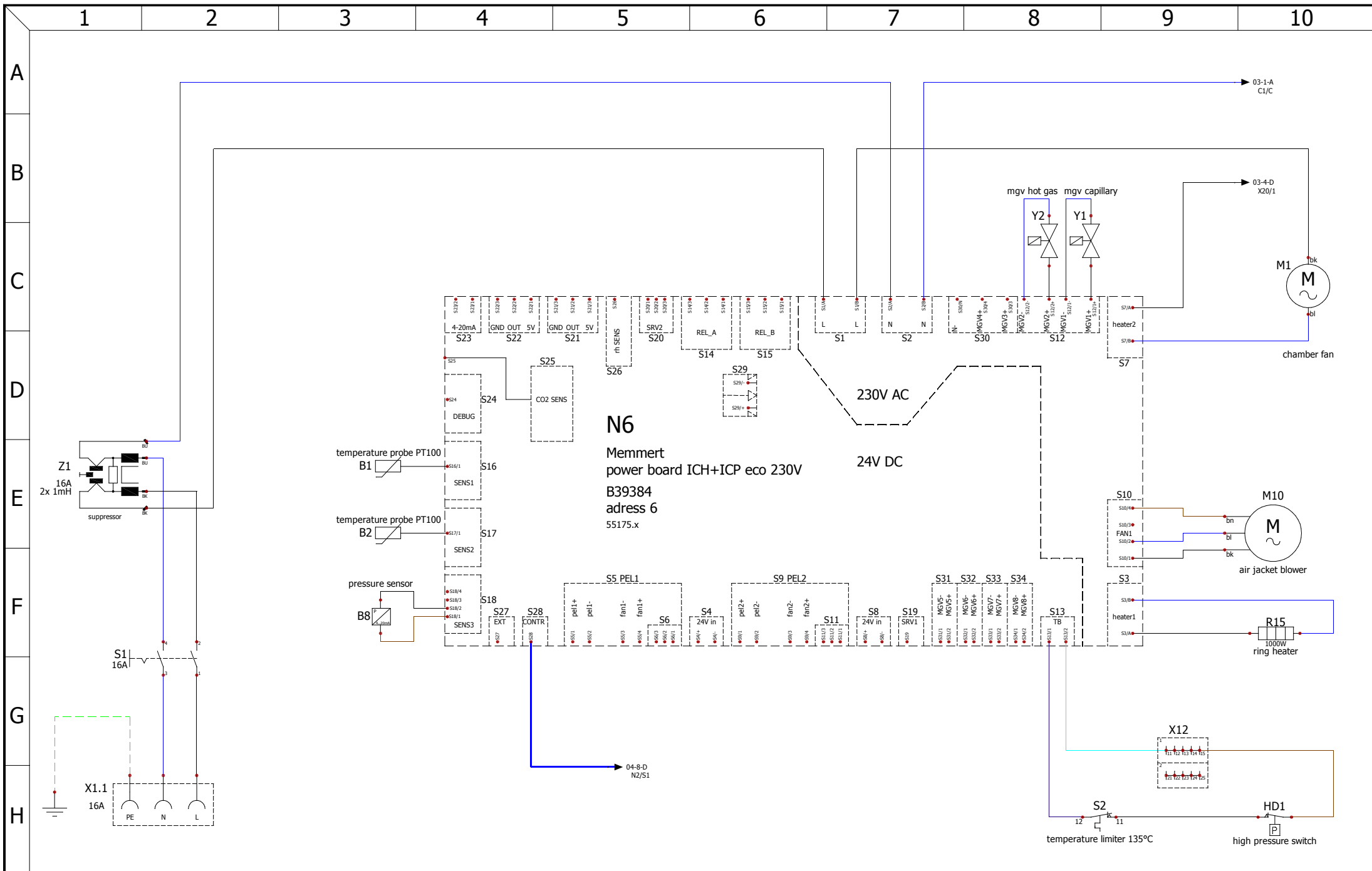


power board 1			designation:	unit or size:	rev	modification	date	name	page
			ICP eco	260					02
created by	date	name	drawing number:	nominal voltage:					08
checked by	31.08.2018	M.Hammerbacher	D39571	230V	A	suppressor Z1 added	07.11.2018	Chr.Seyfarth	

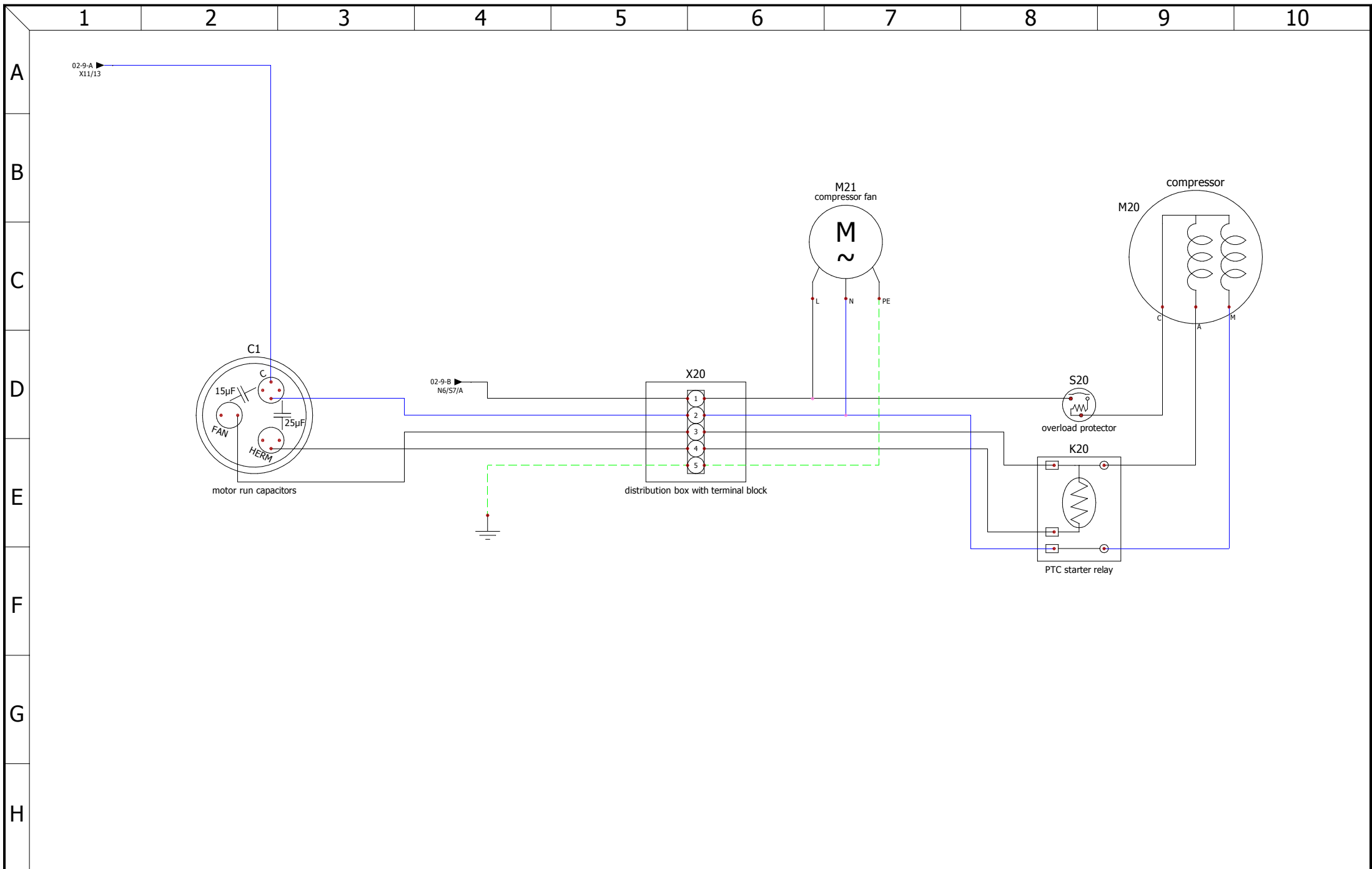



compressor			designation:	unit or size:	rev	modification	date	name	page
			ICP eco	260					03
created by	27.08.2018	Chr.Seyfarth	drawing number:	nominal voltage:					08
checked by	31.08.2018	M.Hammerbacher	D39571	230V	A	suppressor Z1 added	07.11.2018	Chr.Seyfarth	

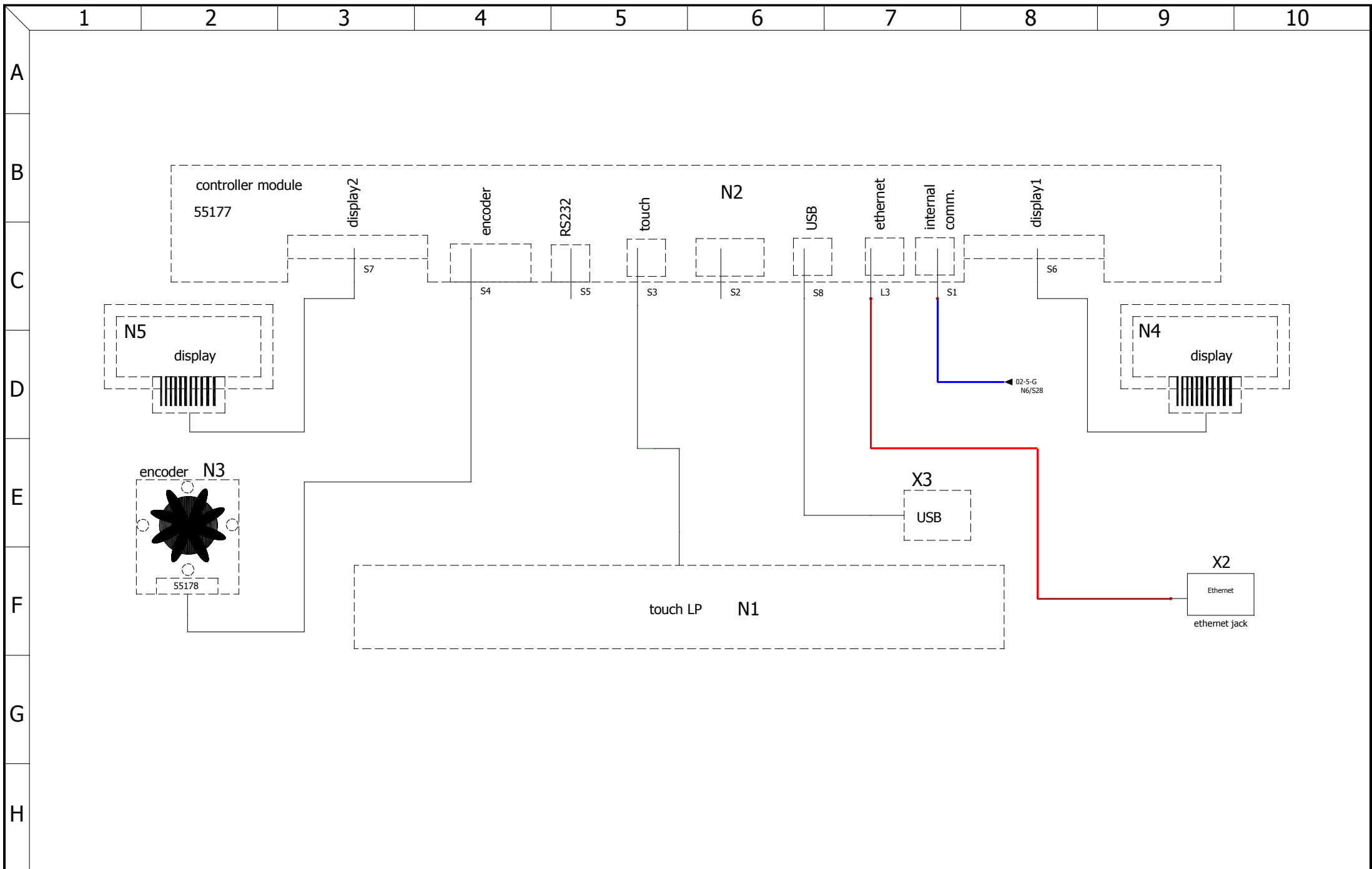

 Experts in Thermostatics	controller unit		designation:	unit or size:	rev	modification	date	name	page
			ICP eco	260					04
	created by	date	name	drawing number:	nominal voltage:				08
	checked by	31.08.2018	M.Hammerbacher	D39571	230V	A	suppressor Z1 added	07.11.2018	Chr.Seyfarth

- hot gas pipe
- liquid gas pipe
- suction pipe
- insulated pipe

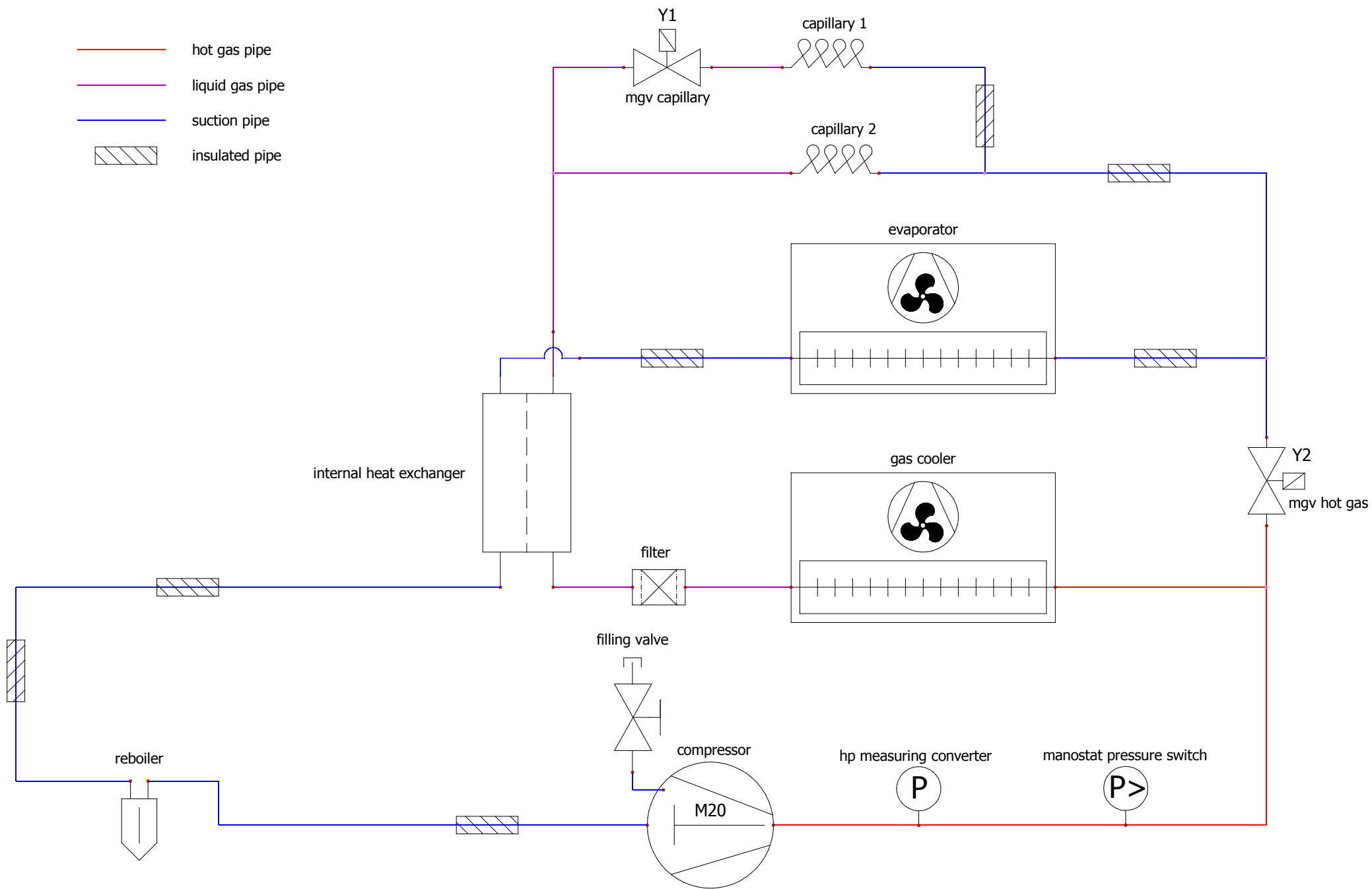

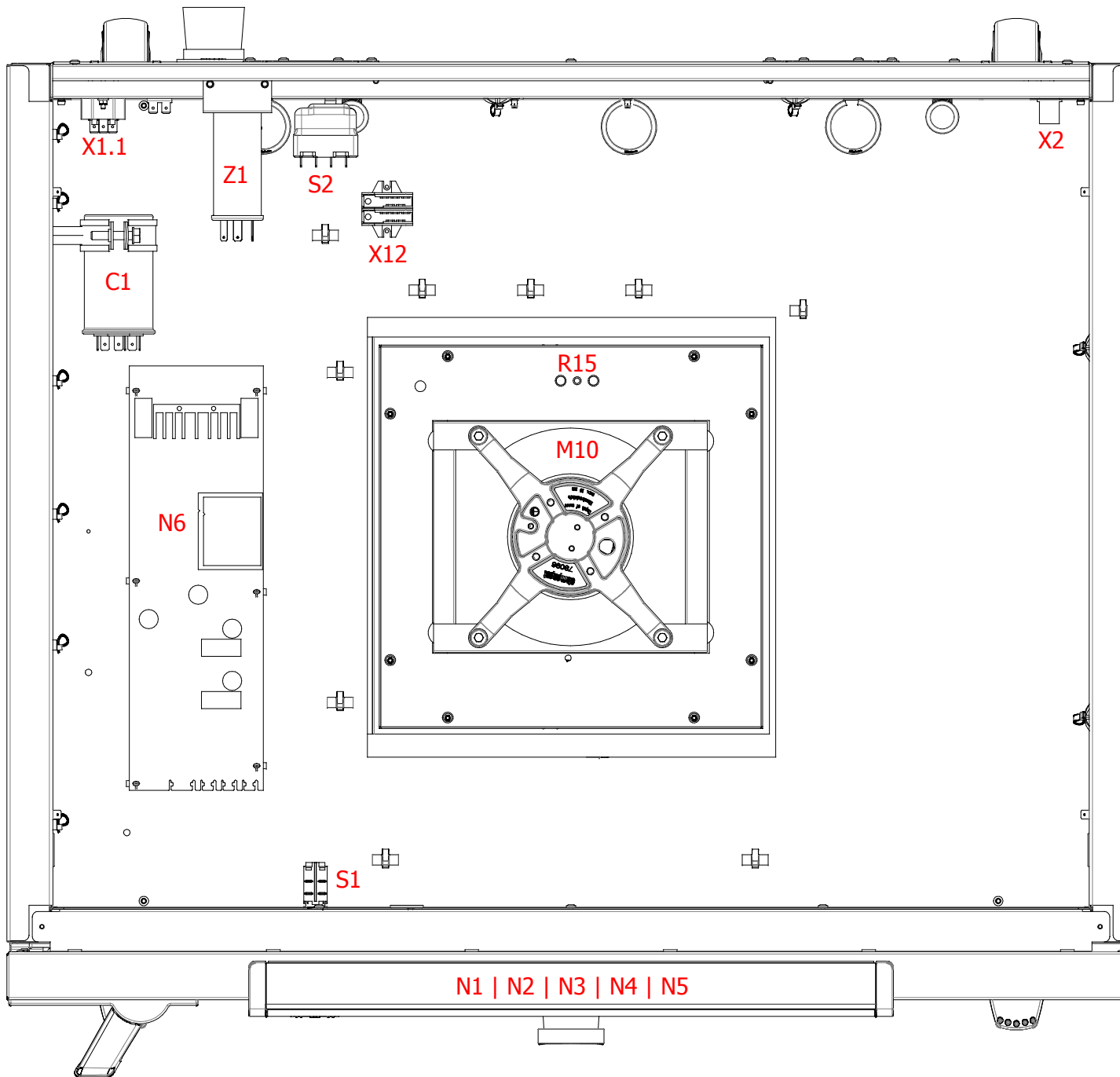


control enclosure		
	date	name
created by	27.08.2018	Chr.Seyfarth
checked by	31.08.2018	M.Hammerbacher

designation:	ICP eco
drawing number:	D39571

unit or size:	260
nominal voltage:	230V

rev	modification	date	name	page
				06
A	suppressor Z1 added	07.11.2018	Chr.Seyfarth	08

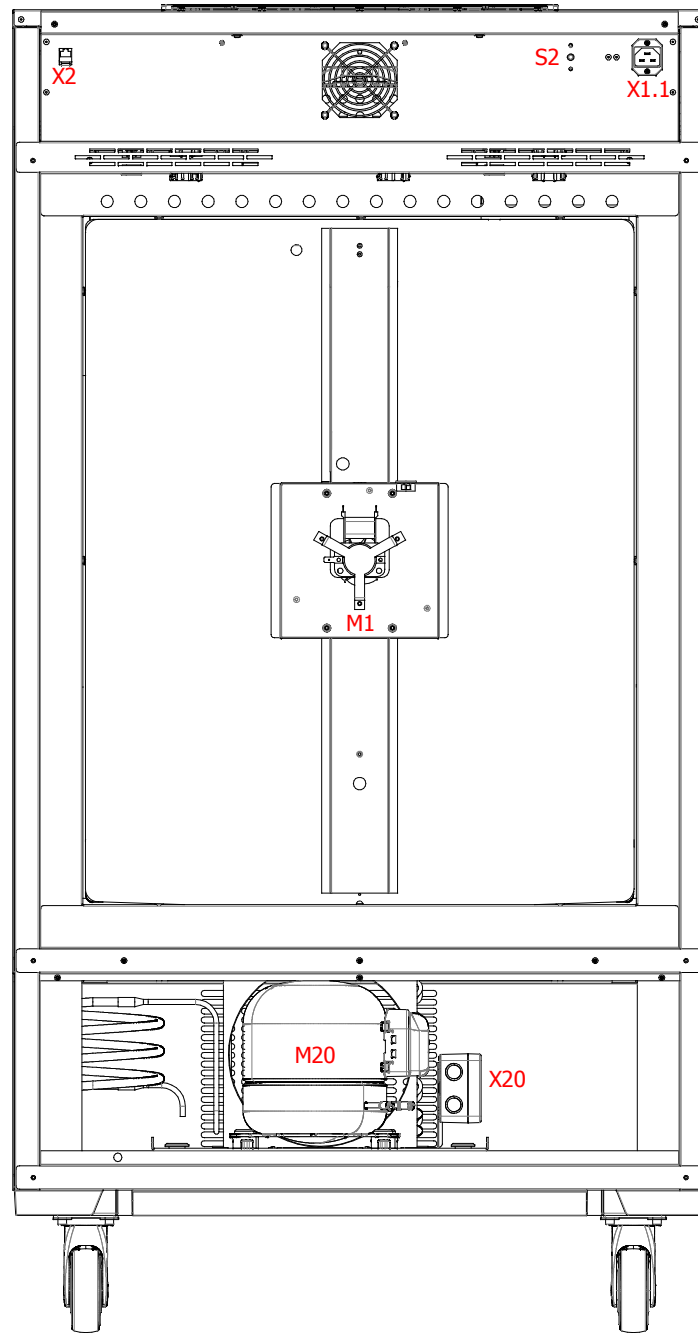

back

	date	name
created by	27.08.2018	Chr.Seyfarth
checked by	31.08.2018	M.Hammerbacher

designation:
ICP eco
drawing number:
D39571

unit or size:
260
nominal voltage:
230V

rev	modification	date	name	page
				07
A	suppressor Z1 added	07.11.2018	Chr.Seyfarth	08

mark	article number	description	quantity
-	B39570	cable harness ICP eco 260	1
-	E07480	compressor unit R744	1
B1 B2	E05952	temperature sensors	2
C1	E07567	motor run capacitors	1
M1	B32012	fan R2K150	1
M10	B24388	fan air jacket ICP800+ ring heating + temperature limiter	1
N1 N2 N3 N4 N5	B30890	controller unit small plus twin display	1
N6	B39384	power board ICH+ICP eco 230V	1
R15	E00119	ring heater	1
S1	E06137	power push-button-switch	1
S2	E02340	temperature limiter 135°C	1
X2	E06251	ethernet jack	1
X12	E07368	distribution block 2pol.	1
X1.1	E06135	connector plug	1
Z1	E07420	interference suppression filter	1